


Prostaglandin E Metabolite EIA Kit

Item No. 514531

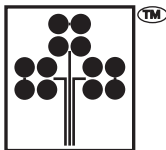

ACE

TABLE OF CONTENTS

GENERAL INFORMATION	3 Materials Supplied
	4 Precautions
	4 If You Have Problems
	4 Storage and Stability
	5 Materials Needed but Not Supplied
INTRODUCTION	6 Background
	6 About This Assay
	8 Description of ACE™ Competitive EIAs
	9 Biochemistry of Acetylcholinesterase
	11 Definition of Key Terms
PRE-ASSAY PREPARATION	12 Buffer Preparation
	14 Sample Preparation
	15 Sample Purification
ASSAY PROTOCOL	16 Derivatization of Standards and Samples to PGEM
	19 Preparation of Assay-Specific Reagents
	21 Plate Set Up
	22 Performing the Assay
ANALYSIS	25 Calculations
	27 Performance Characteristics
RESOURCES	31 Troubleshooting
	32 Additional Reading
	32 References
	33 Related Products
	34 Warranty and Limitation of Remedy
	35 Plate Template
	36 Notes

GENERAL INFORMATION

Materials Supplied

Item Number	Item	96 wells Quantity/Size	480 wells Quantity/Size
414532	Prostaglandin E Metabolite EIA Antiserum	1 vial/100 dtn	1 vial/500 dtn
414530	Prostaglandin E Metabolite AChE Tracer	1 vial/100 dtn	1 vial/500 dtn
414534	Prostaglandin E Metabolite EIA Standard	1 vial	1 vial
400060	EIA Buffer Concentrate (10X)	2 vials/10 ml	4 vials/10 ml
400062	Wash Buffer Concentrate (400X)	1 vial/5 ml	1 vial/12.5 ml
400035	Polysorbate 20	1 vial/3 ml	1 vial/3 ml
400005/400007	Mouse Anti-Rabbit IgG Coated Plate	1 plate	5 plates
400012	96-Well Cover Sheet	1 cover	5 covers
400050	Ellman's Reagent	3 vials/100 dtn	6 vials/250 dtn
400032	Phosphate Buffer	1 vial/100 dtn	2 vials/250 dtn
400027	Carbonate Buffer	1 vial/100 dtn	1 vial/500 dtn
400040	EIA Tracer Dye	1 vial	1 vial
400042	EIA Antiserum Dye	1 vial	1 vial

If any of the items listed above are damaged or missing, please contact our Customer Service department at (800) 364-9897 or (734) 975-3999. We cannot accept any returns without prior authorization.



WARNING: This product is for laboratory research use only: not for administration to humans. Not for human or veterinary diagnostic or therapeutic use.

Precautions

Please read these instructions carefully before beginning this assay.

The reagents in this kit have been tested and formulated to work exclusively with Cayman Chemical's ACE™ EIA Kits. This kit may not perform as described if any reagent or procedure is replaced or modified.

For research use only. Not for human or diagnostic use.

If You Have Problems

Technical Service Contact Information

Phone: 888-526-5351 (USA and Canada only) or 734-975-3888

Fax: 734-971-3641

Email: techserv@caymanchem.com

Hours: M-F 8:00 AM to 5:30 PM EST

In order for our staff to assist you quickly and efficiently, please be ready to supply the lot number of the kit (found on the outside of the box).

Storage and Stability

This kit will perform as specified if stored as directed at -20°C and used before the expiration date indicated on the outside of the box.

Materials Needed But Not Supplied

1. A plate reader capable of measuring absorbance between 405-420 nm.
2. Adjustable pipettes and a repeat pipettor.
3. A source of 'UltraPure' water. Water used to prepare all EIA reagents and buffers must be deionized and free of trace organic contaminants ('UltraPure'). Use activated carbon filter cartridges or other organic scavengers. Glass distilled water (even if double distilled), HPLC-grade water, and sterile water (for injections) are not adequate for EIA. *NOTE: UltraPure water is available for purchase from Cayman (Item No. 400000).*
4. Materials used for **Sample Preparation** (see page 14).

Background

Prostaglandin E₂ (PGE₂) is produced by a variety of cell types which, in general, do not contain the enzymes required for metabolism of PGE₂. Thus, cultured endothelial cells or osteoblasts will release PGE₂ into the culture medium where it will accumulate without appreciable metabolism. The direct assay of PGE₂ from the medium is a good way to measure PGE₂ production from these cells.

PGE₂ is rapidly converted *in vivo* to its 13,14-dihydro-15-keto metabolite (see Figure 1, on page 7), with more than 90% of circulating PGE₂ cleared by a single passage through the lungs. Unfortunately, this metabolite is not chemically stable and undergoes a variable amount of degradation to PGA products. For this reason, blood, urine, or other samples from whole animals or humans often contain very little intact PGE₂, and measurement of the metabolites is necessary to provide a reliable estimate of actual PGE₂ production.

About This Assay

Cayman's Prostaglandin E Metabolite (PGEM) assay is a competitive assay that converts 13,14-dihydro-15-keto PGA₂ and 13,14-dihydro-15-keto PGE₂ to a single, stable derivative that can be easily quantified. This assay is, therefore, the method of choice if the samples in question have undergone extensive metabolism prior to collection. The EIA typically displays an IC₅₀ (50% B/B₀) value of approximately 10 pg/ml and a detection limit (80% B/B₀) of approximately 2 pg/ml.

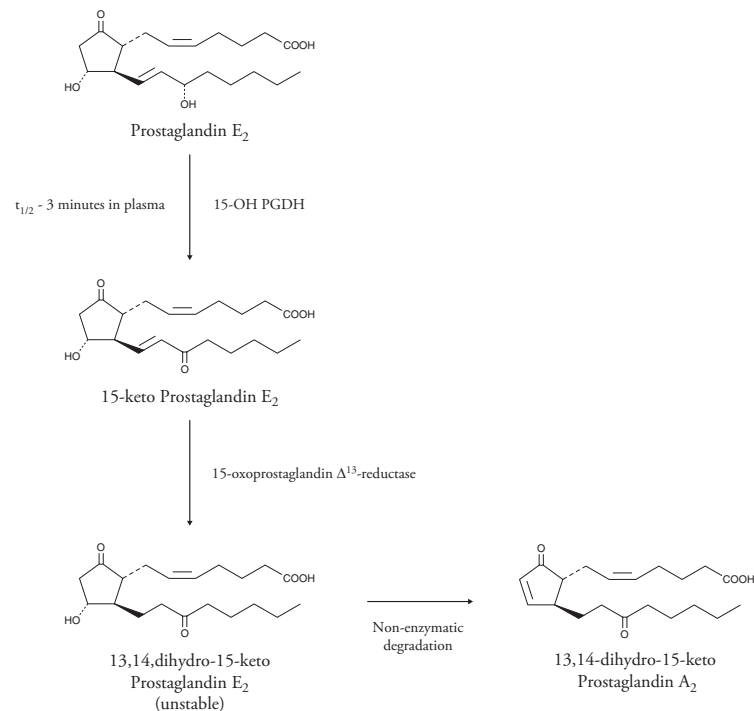


Figure 1. Metabolism of PGE₂

Description of ACE™ Competitive EIAs^{1,2}

This assay is based on the competition between Prostaglandin E Metabolite (PGEM) and a PGEM-acetylcholinesterase (AChE) conjugate (PGEM Tracer) for a limited number of PGEM-specific rabbit antiserum binding sites. Because the concentration of the PGEM Tracer is held constant while the concentration of PGEM varies, the amount of PGEM Tracer that is able to bind to the rabbit antiserum will be inversely proportional to the concentration of PGEM in the well. This rabbit antiserum-PGEM (either free or tracer) complex binds to the mouse monoclonal anti-rabbit IgG that has been previously attached to the well. The plate is washed to remove any unbound reagents, and then Ellman's Reagent (which contains the substrate to AChE) is added to the well. The product of this enzymatic reaction has a distinct yellow color and absorbs strongly at 412 nm. The intensity of this color, determined spectrophotometrically, is proportional to the amount of PGEM Tracer bound to the well, which is inversely proportional to the amount of free PGEM present in the well during the incubation; or

$$\text{Absorbance} \propto [\text{Bound PGEM Tracer}] \propto 1/[\text{PGEM}]$$

A schematic of this process is shown in Figure 2, below.

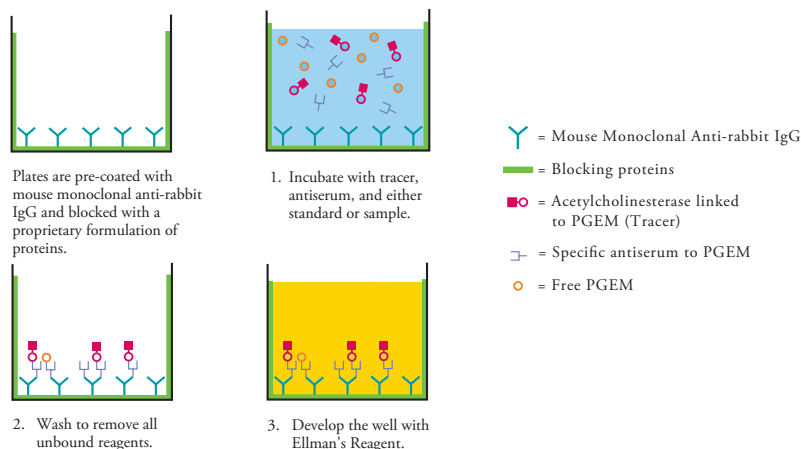


Figure 2. Schematic of the ACE™ EIA

Biochemistry of Acetylcholinesterase

The electric organ of the electric eel, *Electrophorus electricus*, contains an avid acetylcholinesterase (AChE) capable of massive catalytic turnover during the generation of its electrochemical discharges. The electric eel AChE has a clover leaf-shaped tertiary structure consisting of a triad of tetramers attached to a collagen-like structural fibril. This stable enzyme is capable of high turnover ($64,000 \text{ s}^{-1}$) for the hydrolysis of acetylthiocholine.

A molecule of the analyte covalently attached to a molecule of AChE serves as the tracer in ACE™ enzyme immunoassays. Quantification of the tracer is achieved by measuring its AChE activity with Ellman's Reagent. This reagent consists of acetylthiocholine and 5,5'-dithio-*bis*-(2-nitrobenzoic acid). Hydrolysis of acetylthiocholine by AChE produces thiocholine (see Figure 3, on page 10). The non-enzymatic reaction of thiocholine with 5,5'-dithio-*bis*-(2-nitrobenzoic acid) produces 5-thio-2-nitrobenzoic acid, which has a strong absorbance at 412 nm ($\epsilon = 13,600$).

AChE has several advantages over other enzymes commonly used for enzyme immunoassays. Unlike horseradish peroxidase, AChE does not self-inactivate during turnover. This property of AChE also allows re-development of the assay if it is accidentally splashed or spilled. In addition, the enzyme is highly stable under the assay conditions, has a wide pH range (pH 5-10), and is not inhibited by common buffer salts or preservatives. Since AChE is stable during the development step, it is unnecessary to use a 'stop' reagent, and the plate may be read whenever it is convenient.

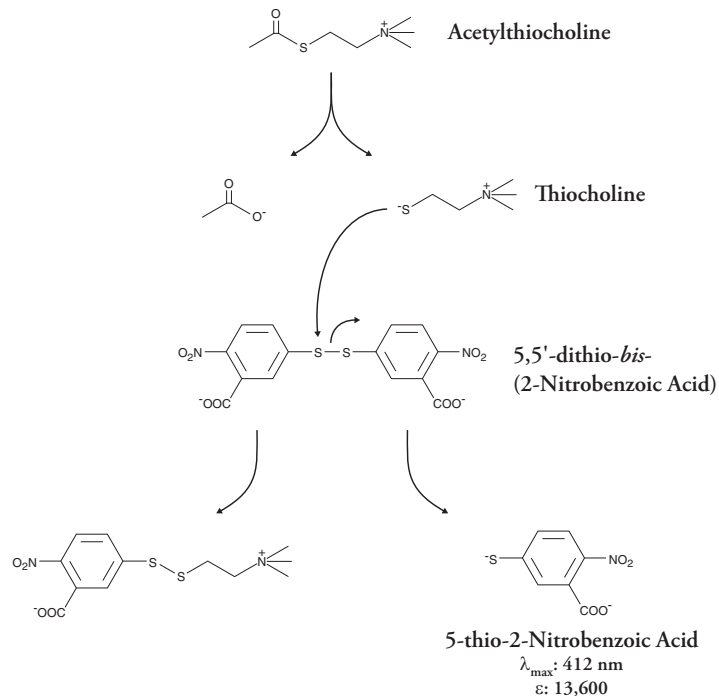


Figure 3. Reaction catalyzed by acetylcholinesterase

Definition of Key Terms

Blank: background absorbance caused by Ellman's Reagent. The blank absorbance should be subtracted from the absorbance readings of all the other wells.

Total Activity: total enzymatic activity of the AChE-linked tracer. This is analogous to the specific activity of a radioactive tracer.

NSB (Non-Specific Binding): non-immunological binding of the tracer to the well. Even in the absence of specific antibody a very small amount of tracer still binds to the well; the NSB is a measure of this low binding.

B_0 (Maximum Binding): maximum amount of the tracer that the antibody can bind in the absence of free analyte.

%B/ B_0 (%Bound/Maximum Bound): ratio of the absorbance of a particular sample or standard well to that of the maximum binding (B_0) well.

Standard Curve: a plot of the %B/ B_0 values *versus* concentration of a series of wells containing various known amounts of analyte.

Dtn: determination, where one dtn is the amount of reagent used per well.

PRE-ASSAY PREPARATION

NOTE: Water used to prepare all EIA reagents and buffers must be deionized and free of trace organic contaminants ('UltraPure'). Use activated carbon filter cartridges or other organic scavengers. Glass distilled water (even if double distilled), HPLC-grade water, and sterile water (for injections) are not adequate for EIA. UltraPure water may be purchased from Cayman (Item No. 400000).

Buffer Preparation

Store all diluted buffers at 4°C; they will be stable for about two months

1. EIA Buffer Preparation

Dilute the contents of one vial of EIA Buffer Concentrate (10X) (Item No. 400060) with 90 ml of UltraPure water. Be certain to rinse the vial to remove any salts that may have precipitated. *NOTE: It is normal for the concentrated buffer to contain crystalline salts after thawing. These will completely dissolve upon dilution with water.*

2. Wash Buffer Preparation

5 ml vial Wash Buffer Concentrate (400X) (96-well kit; Item No. 400062):
Dilute to a total volume of 2 liters with UltraPure water and add 1 ml of Polysorbate 20 (Item No. 400035).

OR

12.5 ml vial Wash Buffer Concentrate (400X) (480-well kit; Item No. 400062):
Dilute to a total volume of 5 liters with UltraPure water and add 2.5 ml of Polysorbate 20 (Item No. 400035).

Smaller volumes of Wash Buffer can be prepared by diluting the Wash Buffer Concentrate 1:400 and adding Polysorbate 20 (0.5 ml/liter of Wash Buffer).

NOTE: Polysorbate 20 is a viscous liquid and cannot be measured by a regular pipette. A positive displacement pipette or a syringe should be used to deliver small quantities accurately.

3. Phosphate Buffer

Prepare a 1 M Phosphate Buffer solution by dissolving the contents of the 100 dtn vial of Phosphate Buffer (Item No. 400032) in 30 ml UltraPure water, or dissolve the contents of one of the 250 dtn vials of Phosphate Buffer (Item No. 400032) in 75 ml UltraPure water.

4. Carbonate Buffer

Prepare a 1 M Carbonate Buffer solution by dissolving the contents of the 100 dtn vial of Carbonate Buffer (Item No. 400027) in 25 ml UltraPure water, or dissolve the contents of the 500 dtn vial of Carbonate Buffer (Item No. 400027) in 125 ml UltraPure water.

5. PGEM Assay Buffer

Prepare 20 ml of PGEM Assay Buffer by combining 13 ml EIA Buffer, 3 ml Carbonate Buffer, and 4 ml Phosphate Buffer. This quantity of buffer should be more than sufficient to complete one 96-well plate.

Sample Preparation

This assay has been validated for a wide range of samples including urine, tissue culture supernatants and plasma. Proper sample storage and preparation are essential for consistent and accurate results. Please read this section carefully before beginning the assay.

General Precautions

- All samples must be free of organic solvents prior to assay.
- Samples should be assayed immediately after collection; samples that cannot be assayed immediately should be stored at -80°C .
- Samples of rabbit origin may contain antibodies which interfere with the assay by binding to the mouse anti-rabbit plate. We recommend that all rabbit samples be purified prior to use in this assay.

Testing for Interference

In general, tissue culture supernatant and urine samples may be derivatized, diluted with PGE Metabolite Buffer and added directly to the assay well. Plasma, serum, and other heterogeneous mixtures such as lavage fluids and aspirates often contain contaminants which can interfere in the assay. It is best to check for interference before embarking on a large number of sample measurements. To test for interference, derivatize one or two test samples, then dilute with PGE Metabolite Buffer to obtain at least two different dilutions of each sample between ~ 2 and 50 pg/ml (*i.e.*, between 20-80% B/B₀). If the two different dilutions of the sample show good correlation (differ by 20% or less) in the final calculated PGEM concentrations, purification is not required. If you do not see good correlation of the different dilutions, purification is advised. The protocol described on page 15 is one such method.

Plasma

Analysis of plasma samples without purification may lead to inconsistent results. We recommend using the purification protocol described below for all plasma samples.

Plasma samples should be collected in vacutainers containing sodium citrate, heparin, or EDTA. Vacutainers can also be supplemented with indomethacin to give a final concentration of at least $10\text{ }\mu\text{M}$. Indomethacin will prevent *ex vivo* formation of prostaglandins, which have the potential to interfere with this assay (although most prostaglandins do not appear to exhibit any cross-reactivity (see page 30)). Samples that cannot be assayed immediately should be stored at -80°C .

Sample Purification

Acetone Precipitation

Acetone precipitation is recommended for removal of proteins from heterogeneous samples. For some samples, acetone precipitation may be the only purification that is required.

1. Aliquot samples into clean tubes.
2. Add four volumes of ice-cold acetone.
3. Incubate at -20°C for a minimum of thirty minutes.
4. Pellet proteins by centrifugation in a tabletop centrifuge at a minimum of 400 x g for five minutes.
5. Transfer supernatants to clean tubes.
6. Dry under a gentle stream of nitrogen.
7. Resuspend samples in the original volume of EIA Buffer.
8. The samples are now ready for derivatization

Derivatization of Standards and Samples to PGEM

Derivatization Hints

- Allow the derivatization to proceed overnight to ensure that all the PGE₂ metabolites derivatize completely.
- Derivatize all standards and samples for the same amount of time.

Derivatization of the PGEM EIA Standard

Equilibrate a pipette tip in ethanol by repeatedly filling and expelling the tip with ethanol several times. Using the equilibrated pipette tip, transfer 100 µl of the PGEM EIA Standard (Item No. 414534) into a clean test tube, then dilute with 900 µl UltraPure water. The concentration of this solution (the bulk standard) will be 40 ng/ml.

Aliquot 50 µl of this solution into a clean tube and dilute to a total volume of 1 ml with EIA Buffer (*i.e.*, add 950 µl). Add 300 µl of Carbonate Buffer and incubate at 37°C overnight. Then add 400 µl Phosphate Buffer and 300 µl EIA Buffer. This solution is 1,000 pg/ml.

Derivatization of the Samples

Aliquot 500 µl of each sample into a clean test tube. Add 150 µl of Carbonate Buffer and incubate overnight at 37°C. Then add 200 µl of Phosphate Buffer and 150 µl of EIA Buffer. Depending on your sample type, the samples may now be ready to use directly in the assay.

NOTE: Because of the high salt concentration in the PGEM Standards, it is crucial that PGEM Assay Buffer is used rather than EIA Buffer to dilute samples if further dilution is necessary.

For human plasma samples, the following ethyl acetate extraction protocol is suggested after derivatization.

Acidification and Ethyl Acetate Extraction

We suggest the following procedure for samples requiring additional purification following derivatization:

1. Acidify the sample to ~pH 4 by the addition of 1M acetate buffer. (To avoid having to measure the pH of each individual sample, adjust the pH of an equivalent volume of sample matrix to pH 4.0 using the 1M acetate buffer. *NOTE: For samples of different volumes, the amount of acid should be adjusted to maintain this ratio of acid to sample.*)
2. Add two volumes of ethyl acetate to each derivatized sample.
3. Vortex thoroughly to mix.
4. Centrifuge samples at a minimum of 200 x g for five minutes
5. Transfer the ethyl acetate (upper layer) to a clean glass tube.
6. Repeat this extraction three more times, combining all of the ethyl acetate layers from each individual sample.
7. Dry samples under a gentle stream of nitrogen.
8. Resuspend samples in PGEM Assay Buffer.
9. The samples are now ready to use in the assay

NOTE: Because of the high salt concentration in the PGEM Standards, it is crucial that PGEM Assay Buffer is used rather than EIA Buffer to dilute samples if further dilution is necessary.

Preparing the Standard Curve

NOTE: Because of the high salt concentration in the 1,000 pg/ml solution, all the points of the standard curve must contain the same salt concentration. Thus, when performing the serial dilution, use the PGEM Assay Buffer.

To prepare the standard for use in EIA: Obtain eight clean test tubes and number them #1 through #8. Aliquot 950 μ l PGEM Assay Buffer to tube #1 and 500 μ l PGEM Assay Buffer to tubes #2-8. Transfer 50 μ l of the derivatized standard (1,000 pg/ml) to tube #1 and mix thoroughly. Serially dilute the standard by removing 500 μ l from tube #1 and placing it into tube #2; mix thoroughly. Next, remove 500 μ l from tube #2 and place it into tube #3; mix thoroughly. Repeat this process for tubes #4-8. These diluted standards should not be stored for more than 24 hours.

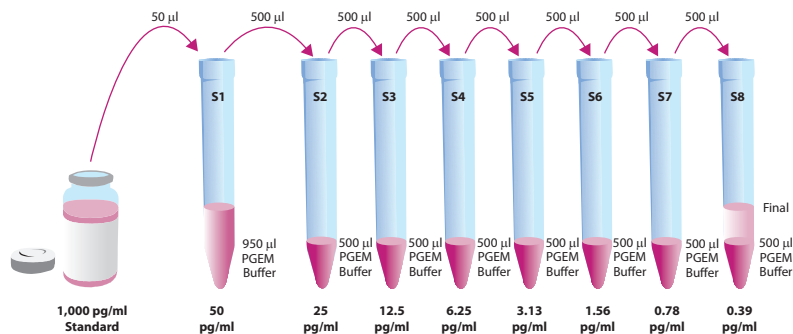


Figure 4. Preparation of the PGEM standards

Preparation of Assay-Specific Reagents

PGEM AChE Tracer

Reconstitute the PGEM AChE Tracer as follows:

100 dtn PGEM AChE Tracer (96-well kit; Item No. 414530): Reconstitute with 6 ml EIA Buffer.

OR

500 dtn PGEM AChE Tracer (480-well kit; Item No. 414530): Reconstitute with 30 ml EIA Buffer.

Store the reconstituted PGEM AChE Tracer at 4°C (*do not freeze!*) and use within four weeks. A 20% surplus of tracer has been included to account for any incidental losses.

Tracer Dye Instructions (optional)

This dye may be added to the tracer, if desired, to aid in visualization of tracer-containing wells. Add the dye to the reconstituted tracer at a final dilution of 1:100 (add 60 μ l of dye to 6 ml tracer or add 300 μ l of dye to 30 ml of tracer).

NOTE: Do not store tracer with dye for more than 24 hours.

PGEM EIA Antiserum

Reconstitute the PGEM EIA Antiserum as follows:

100 dtn PGEM EIA Antiserum (96-well kit; Item No. 414532): Reconstitute with 6 ml EIA Buffer.

OR

500 dtn PGEM EIA Antiserum (480-well kit; Item No. 414532): Reconstitute with 30 ml EIA Buffer.

Store the reconstituted PGEM EIA Antiserum at 4°C. It will be stable for at least four weeks. A 20% surplus of antiserum has been included to account for any incidental losses.

Antiserum Dye Instructions (optional)

This dye may be added to the antiserum, if desired, to aid in visualization of antiserum-containing wells. Add the dye to the reconstituted antiserum at a final dilution of 1:100 (add 60 µl of dye to 6 ml antiserum or add 300 µl of dye to 30 ml of antiserum). *NOTE: Do not store antiserum with dye for more than 24 hours.*

Plate Set Up

The 96-well plate(s) included with this kit is supplied ready to use. It is not necessary to rinse the plate(s) prior to adding the reagents. *NOTE: If you do not need to use all the strips at once, place the unused strips back in the plate packet and store at 4°C. Be sure the packet is sealed with the desiccant inside.*

Each plate or set of strips must contain a minimum of two blanks (Blk), two non-specific binding wells (NSB), two maximum binding wells (B_0), and an eight point standard curve run in duplicate. *NOTE: Each assay must contain this minimum configuration in order to ensure accurate and reproducible results.* Each sample should be assayed at two dilutions and each dilution should be assayed in duplicate. For statistical purposes, we recommend assaying samples in triplicate.

A suggested plate format is shown in Figure 5, below. The user may vary the location and type of wells present as necessary for each particular experiment. The plate format provided below has been designed to allow for easy data analysis using a convenient spreadsheet offered by Cayman (see page 26, for more details). We suggest you record the contents of each well on the template sheet provided (see page 35).

	1	2	3	4	5	6	7	8	9	10	11	12
A	Blk	S1	S1	1	1	1	9	9	9	17	17	17
B	Blk	S2	S2	2	2	2	10	10	10	18	18	18
C	NSB	S3	S3	3	3	3	11	11	11	19	19	19
D	NSB	S4	S4	4	4	4	12	12	12	20	20	20
E	B_0	S5	S5	5	5	5	13	13	13	21	21	21
F	B_0	S6	S6	6	6	6	14	14	14	22	22	22
G	B_0	S7	S7	7	7	7	15	15	15	23	23	23
H	TA	S8	S8	8	8	8	16	16	16	24	24	24

Blk - Blank
TA - Total Activity
NSB - Non-Specific Binding
 B_0 - Maximum Binding
S1-S8 - Standards 1-8
1-24 - Samples

Figure 5. Sample plate format

Performing the Assay

Pipetting Hints

- Use different tips to pipette each reagent.
- Before pipetting each reagent, equilibrate the pipette tip in that reagent (*i.e.*, slowly fill the tip and gently expel the contents, repeat several times).
- Do not expose the pipette tip to the reagent(s) already in the well.

Addition of the Reagents

1. PGEM Buffer

Add 50 μ l EIA Buffer and 50 μ l of PGEM Buffer to Non-Specific Binding (NSB) wells. Add 50 μ l PGEM Buffer to Maximum Binding (B_0) wells.

2. PGEM EIA Standard

Add 50 μ l from tube #8 to both of the lowest standard wells (S8). Add 50 μ l from tube #7 to each of the next two standard wells (S7). Continue with this procedure until all the standards are aliquoted. The same pipette tip should be used to aliquot all the standards. Before pipetting each standard, be sure to equilibrate the pipette tip in that standard.

3. Samples

Add 50 μ l of sample per well. Each sample should be assayed at a minimum of two dilutions. Each dilution should be assayed in duplicate (triplicate recommended).

4. PGEM AChE Tracer

Add 50 μ l to each well *except* the Total Activity (TA) and the Blank (Blk) wells.

5. PGEM EIA Antiserum

Add 50 μ l to each well *except* the Total Activity (TA), the Non-Specific Binding (NSB), and the Blank (Blk) wells.

Well	PGEM Buffer	EIA Buffer	Standard/ Sample	Tracer	Antibody
Blk	-	-	-	-	-
TA	-	-	-	5 μ l (at devel. step)	-
NSB	50 μ l	50 μ l	-	50 μ l	-
B_0	50 μ l	-	-	50 μ l	50 μ l
Std/Sample	-	-	50 μ l	50 μ l	50 μ l

Table 1. Pipetting summary

Incubation of the Plate

Cover each plate with plastic film (Item No. 400012) and incubate for 18 hours at room temperature.

Development of the Plate

1. Reconstitute Ellman's Reagent immediately before use (20 ml of reagent is sufficient to develop 100 wells):

100 dtn vial Ellman's Reagent (96-well kit; Item No. 400050): Reconstitute with 20 ml of UltraPure water.

OR

250 dtn vial Ellman's Reagent (480-well kit; Item No. 400050): Reconstitute with 50 ml of UltraPure water.

NOTE: Reconstituted Ellman's Reagent is unstable and should be used the same day it is prepared; protect the Ellman's Reagent from light when not in use. Extra vials of the reagent have been provided should a plate need to be re-developed or multiple assays be run on different days.

2. Empty the wells and rinse five times with Wash Buffer.
3. Add 200 μl of Ellman's Reagent to each well
4. Add 5 μl of tracer to the Total Activity wells.
5. Cover the plate with plastic film. Optimum development is obtained by using an orbital shaker equipped with a large, flat cover to allow the plate(s) to develop in the dark. This assay typically develops (*i.e.*, B_0 wells ≥ 0.3 A.U. (blank subtracted)) in 60-90 minutes.

Reading the Plate

1. Wipe the bottom of the plate with a clean tissue to remove fingerprints, dirt, etc.
2. Remove the plate cover being careful to keep Ellman's Reagent from splashing on the cover. *NOTE: Any loss of Ellman's Reagent will affect the absorbance readings. If Ellman's Reagent is present on the cover, use a pipette to transfer the Ellman's Reagent into the well. If too much Ellman's Reagent has splashed on the cover to easily redistribute back into the wells, wash the plate three times with Wash Buffer and repeat the development with fresh Ellman's Reagent.*
3. Read the plate at a wavelength between 405 and 420 nm. The absorbance may be checked periodically until the B_0 wells have reached a minimum of 0.3 A.U. (blank subtracted). The plate should be read when the absorbance of the B_0 wells in the range of 0.3-1.0 A.U. (blank subtracted). If the absorbance of the wells exceeds 1.5, wash the plate, add fresh Ellman's Reagent and let it develop again.

ANALYSIS

Many plate readers come with data reduction software that plots data automatically. Alternatively a spreadsheet program can be used. The data should be plotted as either $\%B/B_0$ versus log concentration using a four-parameter logistic fit or as logit B/B_0 versus log concentration using a linear fit. *NOTE: Cayman has a computer spreadsheet available for data analysis. Please contact Technical Service or visit our website (www.caymanchem.com/analysis/cia) to obtain a free copy of this convenient data analysis tool.*

Calculations

Preparation of the Data

The following procedure is recommended for preparation of the data prior to graphical analysis.

NOTE: If the plate reader has not subtracted the absorbance readings of the blank wells from the absorbance readings of the rest of the plate, be sure to do that now.

1. Average the absorbance readings from the NSB wells.
2. Average the absorbance readings from the B_0 wells.
3. Subtract the NSB average from the B_0 average. This is the corrected B_0 or corrected maximum binding.
4. Calculate the B/B_0 (Sample or Standard Bound/Maximum Bound) for the remaining wells. To do this, subtract the average NSB absorbance from the S1 absorbance and divide by the corrected B_0 (from Step 3). Repeat for S2-S8 and all sample wells. (To obtain $\%B/B_0$ for a logistic four-parameter fit, multiply these values by 100.)

NOTE: The total activity (TA) values are not used in the standard curve calculations. Rather, they are used as a diagnostic tool; the corrected B_0 divided by the actual TA (10X measured absorbance) will give the % Bound. This value should closely approximate the % Bound that can be calculated from the Sample Data (see page 28). Erratic absorbance values and a low (or no) % Bound could indicate the presence of organic solvents in the buffer or other technical problems (see page 32 for Troubleshooting).

Plot the Standard Curve

Plot %B/B₀ for standards S1-S8 *versus* PGEM concentration using linear (y) and log (x) axes and perform a 4-parameter logistic fit.

Alternative Plot - The data can also be linearized using a logit transformation. The equation for this conversion is shown below. *NOTE: Do not use %B/B₀ in this calculation.*

$$\text{logit} (B/B_0) = \ln [B/B_0 / (1 - B/B_0)]$$

Plot the data as logit (B/B₀) *versus* log concentrations and perform a linear regression fit.

Determine the Sample Concentration

Calculate the B/B₀ (or %B/B₀) value for each sample. Determine the concentration of each sample using the equation obtained from the standard curve. *NOTE: Remember to account for any concentration or dilution of the sample prior to the addition to the well.* Samples with %B/B₀ values greater than 80% or less than 20% should be re-assayed as they generally fall out of the linear range of the standard curve. A 20% or greater disparity between the apparent concentration of two different dilutions of the same sample indicates interference which could be eliminated by purification.

Performance Characteristics

Sample Data

The standard curve presented here is an example of the data typically produced with this kit; however, your results will not be identical to these. You **must** run a new standard curve. Do not use the data below to determine the values of your samples. Your results could differ substantially.

	Raw Data		Average	Corrected
Total Activity	1.752	1.965	1.859	
NSB	-0.001	-0.003	-0.002	
B₀	0.674	0.726		
	0.684	0.719	0.701	0.703

Dose (pg/ml)	Raw Data		Corrected		%B/B ₀	
50	0.138	0.120	0.140	0.122	19.88	17.44
25	0.230	0.213	0.232	0.215	33.08	30.60
12.5	0.335	0.307	0.337	0.309	47.96	44.01
6.25	0.441	0.431	0.443	0.433	63.00	61.63
3.13	0.518	0.521	0.520	0.523	74.00	74.37
1.56	0.592	0.582	0.594	0.584	84.54	83.16
0.78	0.640	0.618	0.642	0.620	91.35	88.20
0.39	0.649	0.687	0.651	0.689	92.58	97.98

Table 2. Typical results

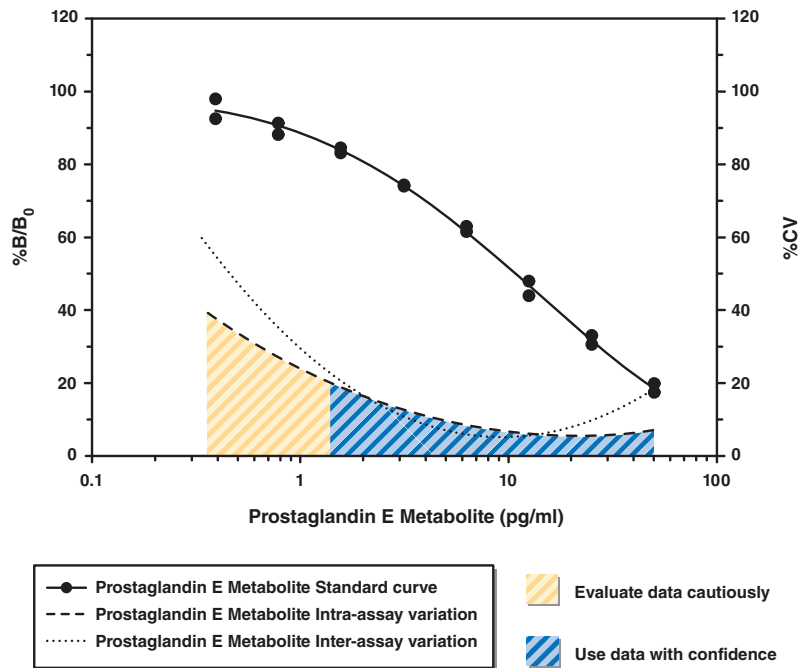


Figure 6. Typical standard curve

Precision:

The intra- and inter-assay CVs have been determined at multiple points on the standard curve. These data are summarized in the graph on page 28 and in the table below.

Dose (pg/ml)	%CV* Intra-assay variation	%CV* Inter-assay variation
50	8.1	18.2
25	5.4	7.2
12.5	5.9	8.3
6.25	5.5	11.2
3.13	12.8	8.0
1.56	25.1	13.4
0.78	23.7	39.3
0.39	N.D.	123

Table 3. Intra- and inter-assay Variation

*%CV represents the variation in concentration (not absorbance) of 40 repetitions of each point on the standard curve as determined using a reference standard curve.

Specificity:

Compound	Cross Reactivity
13,14-dihydro-15-keto PGE ₁ *	100%
13,14-dihydro-15-keto PGE ₂ *	100%
Bicyclo Prostaglandin E ₁	38%
13,14-dihydro-15-keto PGD ₂ *	0.08%
13,14-dihydro-15-keto PGF _{2α} *	0.02%
Arachidonic Acid	<0.01%
Leukotriene B ₄	<0.01%
tetranor-PGEM	<0.01%
tetranor-PGFM	<0.01%
Prostaglandin D ₂	<0.01%
Prostaglandin E ₁	<0.01%
6-keto Prostaglandin E ₁	<0.01%
Prostaglandin E ₂	<0.01%
Prostaglandin F _{1α}	<0.01%
6-keto Prostaglandin F _{1α}	<0.01%
Prostaglandin F _{2α}	<0.01%
Thromboxane B ₂	<0.01%

Table 4. Specificity of the PGEM EIA Antiserum

*Derivatized per Assay Protocol

RESOURCES

Troubleshooting

Problem	Possible Causes	Recommended Solutions
Erratic values; dispersion of duplicates	A. Trace organic contaminants in the water source B. Poor pipetting/technique	A. Replace activated carbon filter or change source of UltraPure water
High NSB (>0.035)	A. Poor washing B. Exposure of NSB wells to specific antibody	A. Re-wash plate and redevelop
Very low B ₀	A. Trace organic contaminants in the water source B. Plate requires additional development time C. Dilution error in preparing reagents	A. Replace activated carbon filter or change source of UltraPure water B. Return plate to shaker and re-read later
Low sensitivity (shift in dose response curve)	Standard is degraded	Replace standard
Analyses of two dilutions of a biological sample do not agree (<i>i.e.</i> , more than 20% difference)	Interfering substances are present	Purify sample prior to analysis by EIA ³
Only Total Activity (TA) wells develop	Trace organic contaminants in the water source	Replace activated carbon filter or change source of UltraPure water

Additional Reading

Go to www.caymanchem.com/514531/references for a list of publications citing the use of Cayman's PGEM Metabolite EIA Kit.

References

1. Maclouf, J., Grassi, J., and Pradelles, P. Development of enzyme-immunoassay techniques for the measurement of eicosanoids, Chapter 5, *in* Prostaglandin and Lipid Metabolism in Radiation Injury. Walden, T.L., Jr. and Hughes, H.N., editors, Plenum Press, Rockville, 355-364 (1987).
2. Pradelles, P., Grassi, J. and Maclouf, J.A. Enzyme immunoassays of eicosanoids using acetylcholinesterase as label: An alternative to radioimmunoassay. *Anal. Chem.* **57**, 1170-1173 (1985).
3. Maxey, K.M., Maddipati, K.R. and Birkmeier, J. Interference in enzyme immunoassays. *J. Clin. Immunoassay* **15**,116-120 (1992).

Related Products

Prostaglandin E₂ EIA Kit - Monoclonal - Item No. 514010
Prostaglandin E₂ Express EIA Kit - Item No. 500141
Prostaglandin Screening EIA Kit - Item No. 514012
SPE Cartridges (C-18) - Item No. 400020
STAT-Prostaglandin E₂ EIA Kit - Item No. 514131
UltraPure Water - Item No. 400000

NOTES

This document is copyrighted. All rights are reserved. This document may not, in whole or part, be copied, photocopied, reproduced, translated, or reduced to any electronic medium or machine-readable form without prior consent, in writing, from Cayman Chemical Company.

©06/03/2011, Cayman Chemical Company, Ann Arbor, MI, All rights reserved. Printed in U.S.A.