

Product Information Sheet

Order: # M1003

Mobicol "Classic"

SUMMARY

shipped at RT; store at RT

Mobicol "Classic" is a 1 ml column with particular handling advantages.



Mobicol "Classic"

- is compatible to the laboratory standard,
- can be centrifuged in microfuges,
- smallest and largest volumes can be treated easily.

Filters of the pore sizes 10 µm, 35 µm or 90 µm are available for Mobicols. The filters can be inserted easily with the help of a plastic tool which is delivered with the columns. The upper large filter fits tightly over the whole length of the column body.

Mobicols can be closed tightly with a screw-on cap and a plug at the outlet. The volume is 800 µl with a small filter and 700 µl with a large filter inserted. Mobicols can be autoclaved at 110 °C for 10 min.

Mobicol "Classic"

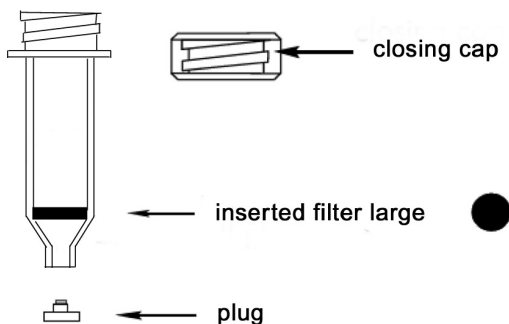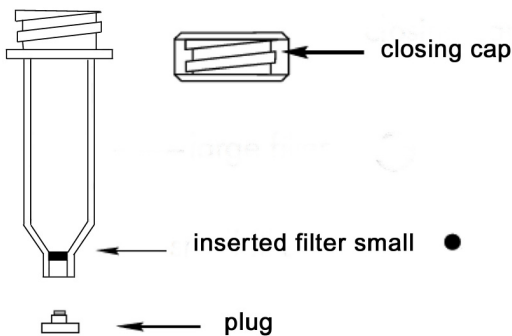


Mobicols can be centrifuged (fit into 1.5 ml or 2 ml tubes)

Product Information Sheet

Order: # M1003

Mobicol "Classic"

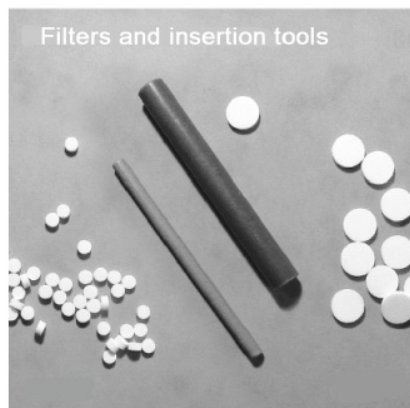

The Mobicol "Classic" M1003 is supplied with a screw cap and a plug for the outlet. Small and large filters with different diameters (10 μm , 35 μm and 90 μm) are separately available, so you can use the filter of your choice. With this features the Mobicol "Classic" can be used for a variety of applications.

Order Information, Shipping and Storage

| Order# | Product | Quantity |
|----------------------------|--|----------|
| M1002 | Mobicol "Classic" with 2 different screw caps, without filters | 50 |
| M1003 | Mobicol "Classic" with 1 closed screw cap and plug, without filters | 50 |
| M2110 | Filter (small) 10 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M513515 | Filter (small) 35 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M2190 | Filter (small) 90 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M2210 | Filter (large) 10 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M523515 | Filter (large) 35 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| M2290 | Filter (large) 90 μm pore size, for Mobicol M1002, M1003, M1050 & M1053 | 50 |
| shipped at RT; store at RT | | |

Contact and Support

MoBiTec GmbH ♦ Lotzestrasse 22a ♦ D-37083 Goettingen ♦ Germany

Customer Service – General inquiries & orders

phone: +49 (0)551 707 22 0
 fax: +49 (0)551 707 22 22
 e-mail: order@mobitec.com

Technical Service – Product information

phone: +49 (0)551 707 22 70
 fax: +49 (0)551 707 22 77
 e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at

www.mobitec.com