

For research use only.

1. About BIOMEDAL LYTAG System.

LYTAG is a new generation of products designed for the expression and purification of proteins in E. coli. The new system is based on a reduced version of the affinity chromatography purification tag C-LYTAG. LYTAG has shorter amino acid sequence (136 aa, 15.21 KDa) than C-LYTAG.

LYTAG, as well as C-LYTAG, specically interacts with LYTRAP resin, a simple, efficient and low cost chromatogreaphic support.

The new LYTAG incorporates an alpha helix polypeptidic sequence which acts as a spacing element between LYTAG and the fused protein, reducing the potencial steric interference between both regions and facilitating elimination of the tag by enterokinase digestion.

2. Description.

PRODUCT	CAT.No.	FORMAT
LYTRAP Resin	RS-3302	30 ml
LYTRAP Resin	RS-3316	250 ml

3. General Protocol

LYTRAP is supplied pre-swollen in 20% ethanol. Pack the required amount of resin (8 ml LYTRAP per litre of culture) in an appropriate column and equilibrate with 5-10 volumes of column equilibration buffer (20 mM sodium phosphate pH 7.0). Be sure the buffer replaces completely the ethanol. For LYTAG protein purification follow the protocol described in "C-LYTAG Purification System" (Biomedal, KT-3246)

NOTES:

- a. During shipping and storage, the resin will settle. We recommend thoroughly resuspending it before pipetting.
- b. The column never can be dry. If this happens, you must resuspend the resin in buffer and repeat the packing.



Binding capacity Bead structure Bead size Recommended flow rate pH stability (<2 h) pH stability (> 2 h) Antimicrobial agent Storage

0.5-3 mg/ml 6% highly cross-linked agarose 45-165 m 1 ml/min 1-14 3-12 Ethanol 20% 4°C for long time periods

Product	Format	Cat.No.
LYTRAP resin	30 ml	RS-3302
LYTRAP resin	250 ml	RS-3316
LYTRAP spin columns	10 units	RS-3319
LYTRAP spin columns	25 units	RS-3320
LYTRAP spin columns	100 units	RS-3321
Purification columns 2.5ml (empty)	5 units	RS-3322
C-LYTAG Purification System	1kit (5 purifications)	KT-3246
C-LYTAG Purification System	1 Kit	KT-3412
LYTAG-Spin Plus Kit	1 Kit	KT-3412

For more information, please visit our WEBSITE or contact us at

Biomedal S.L.

Av. Americo Vespucio, n° 5-E, pl 1°
Parque Científico Tecnológico Cartuja 93
41092 - Sevilla - España/Spain
Tel. +34-954081276 Fax. +34-954081279

www.biomedal.com info@biomedal.com Ed. 3 © 2007 Biomedal, S.L



