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ZF Plate Instruction Manual


The ZF plate is a 96 well microplate that efficiently sorts zebrafish.
This enables smooth imaging of the sorted zebrafish.

1. Contents

Check that the following are included.

- ZF Plate (with transparent cover): 10 plates.
- Instruction Manual: 1

2. Component names

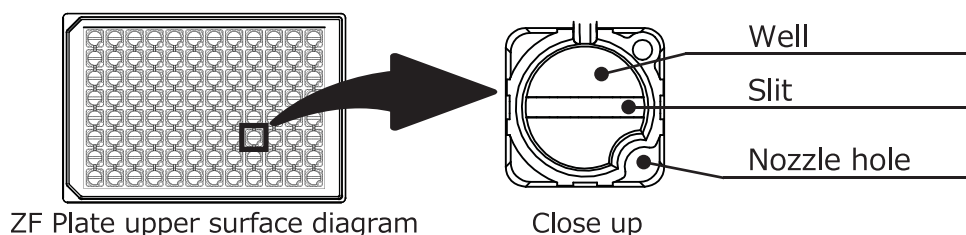3. Procedure

Subject: Anesthetized zebrafish (approx. 3dpf to 6dpf)

Devices used

- ZF Plate
- Micro pipet (100 μ l to 200 μ l and 1 ml, one of each)
- Pipet chip (wide port, low suction type recommended)
- Microplate centrifuge (with flash function)

① Anaesthetized zebrafish are poured into the wells

—Suck 75 μ l (approx.) anesthetic with the zebrafish into the pipet and pour gently into the well.

② Flash processing with centrifuge (spin down)

—Flash process for 1 to 2 turns at 200 to 250Gs of centrifugal force.

③ Taken with imaging device

—After completing imaging, sufficient anesthetic is added to the well and the zebrafish is sucked out with the pipet.

See the URL below for the more details regarding the procedure.

<https://www.hashimoto-inc.co.jp/products/zfプレート/>

4. Specification

Model	HDK-ZFA101-02a	Plate material/color	polystyrene / black
Plate size(mm)	85.48(W) \times 127.76(L) \times 14.35(H)	Bottom material/thickness(mm)	Silica glass / 0.17
Mass(g)	with cover : 72 / without cover : 95	Temperature for use	0 ~ 40 $^{\circ}$ C
Applicable standards	ANSI/SLAS-1 ~ 4 (96well)	Temperature for storage	4 ~ 25 $^{\circ}$ C
Per well maximum capacity(μ l)	250	Sterilized/non-sterilized	non-sterilized

5. Precautions for use

- ⚠ Do not use for applications other than imaging of small fish.
- ⚠ This component is disposable. Do not reuse.
- ⚠ Cultivation of small fish in the well is not possible.
- ⚠ This product cannot be treated with an autoclave.

HASHIMOTO ELECTRONIC INDUSTRY CO.,LTD
3866-12, Takasucho, Matsusaka-shi, Mie, 515-0104, Japan
TEL: +81(0)598-51-3111 FAX : +81(0)598-52-1417
<https://www.hashimoto-inc.co.jp/>