Grid Screen TM **MPD**

User Guide

Grid Screen[™] MPD is a preformulated reagent kit designed to provide a rapid screening method for the crystallization of biological macromolecules. The screen is simple and practical for finding initial crystallization conditions as well as determining the solubility of a macromolecule in MPD between pH 4.0 and 9.0.

Sample Preparation

The macromolecular sample should be homogenous, as pure as is practically possible (>95%) and free of amorphous and particulate material. Remove amorphous material by centrifugation or micro-filtration prior to use (1, 2, 3).

The recommended sample concentration is 5 to 25 mg/ml in water. Initially, the sample should be free of any unnecessary additives in order to observe the effect of the Grid Screen MPD variables. Ideally, the initial screen should be performed with a sample which has been dialyzed against water although ligands, ions, reducing agents, or other additives may be present as required by the sample for solubility, stability, or activity.

Performing the Screen

Since it is the most frequently reported method of crystallization, the following procedure describes the use of Grid Screen MPD with the Hanging Drop Vapor Diffusion method. Grid Screen MPD is also very compatible with the Sitting Drop, Sandwich Drop, Microbatch, and Microdialysis methods. A complete description of the Hanging, Sitting, Sandwich Drop, Dialysis and other crystallization methods are available from the Hampton Research Crystal Growth 101 Library.

1. Prepare a VDX Plate (HR3-140) for Hanging Drop Vapor Diffusion by applying a thin bead of cover slide sealant to the upper edge of each of the 24 reservoirs. One may also use a Greased VDX Plate (HR3-170). Twenty-four reservoirs are to be prepared for a complete Grid Screen MPD. See Figure 1.

Figure 1 Cross section of a reservoir in the VDX plate. Vacuum Grease Well of the VDX Reservoir Crystallization Plate Solution

2. Using a clean pipet tip, pipet 1 ml of Grid Screen MPD reagent A1 into reservoir A1. Discard the pipet tip, add a new pipet tip and pipet 1 ml of Grid Screen MPD reagent A2 into reservoir A2. Repeat the procedure for the remaining 22 Grid Screen MPD reagents using a clean pipet tip for each reagent so as to avoid reagent contamination and carry over.



HR2-215 (pg 1)

3. Pipet 2 µl of the sample to the center of a clean, siliconized 22 mm diameter circle or square cover slide. See Figure 2.

Figure 2



4. Pipet 2 µl of Grid Screen MPD reagent A1 from reservoir A1 into the sample droplet and mix by aspirating and dispensing the droplet several times, keeping the tip in the drop during mixing to avoid foaming. See Figure 2.

5. Working quickly to minimize evaporation, invert the cover slide and droplet over reservoir A1 and seal the cover slide onto the edge of the reservoir. See Figure 3.

Figure 3



6. Repeat operations 3 through 5 for the remaining 23 Grid Screen MPD reagents.

7. If the quantity of sample permits, perform Grid Screen MPD in duplicate and incubate one set of plates at 4°C and the second set at room temperature. Incubate and store the crystallization plates in a stable temperature environment free of vibration.

Examine the Drop

Carefully examine the drops under a stereo microscope (10 to 100x magnification) immediately after setting up the screen. Record all observations and be particularly careful to scan the focal plane for small crystals. Observe the drops once each day for the first week, then once a week there after. Records should indicate whether the drop is clear, contains precipitate, and or crystals. It is helpful to describe the drop contents using descriptive terms. Adding magnitude is also helpful. Example: 4+ yellow/brown fine precipitate, 2+ small bipyramid crystals, clear drop, 3+ needle shaped crystals in 1+ white precipitate. One may also employ a standard numerical scoring scheme (Clear = 0, Precipitate = 1, Crystal = 10, etc). Figure 4 (on page 2) shows typical examples of what one might observe in a crystallization experiment.

Interpreting Grid Screen MPD

Clear drops indicate that either the relative supersaturation of the sample and reagent is too low or the drop has not yet completed equilibration. If

Grid Screen

User Guide

Figure 4

Typical observations in a crystallization experiment



Clear Drop



Skin/Precipitate



Precipitate



Precipitate/Phase



Quasi Crystals



Microcrystals







Rod Cluster



Single Crystal

the drop remains clear after 3 to 4 weeks consider repeating the Grid Screen MPD condition and doubling the sample concentration. If more than 70% Grid Screen MPD drops are clear consider doubling the sample concentration and repeating the entire screen.

Drops containing precipitate indicate that either the relative supersaturation of the sample and reagent is too high, the sample has denatured, or the sample is heterogeneous. To reduce the relative supersaturation, dilute the sample twofold and repeat the Grid Screen MPD condition. If more than 70% Grid Screen MPD drops contain precipitate and no crystals are present, consider diluting the sample concentration in half and repeating the entire screen. If sample denaturation is suspect, take measures to stabilize the sample (add reducing agent, ligands, glycerol, salt, or other stabilizing agents). If the sample is impure, aggregated, or heterogeneous take measures to pursue homogeneity. It is possible to obtain crystals from precipitate so do not discard nor ignore a drop containing precipitate. If possible, examine drops containing precipitate under polarizing optics to differentiate precipitate from microcrystalline material.

If the drop contains a macromolecular crystal the relative supersaturation of the sample and reagent is good. The next step is to optimize the preliminary conditions (pH, salt type, salt concentration, precipitant type, precipitant concentration, sample concentration, temperature, additives, and other crystallization variables) which produced the crystal in order to improve crystal size and quality.

Compare the observations between the 4°C and room temperature incubation to determine the effect of temperature on sample solubility. Different results in the same drops at different temperatures indicate that sample solubility is temperature dependent and that one should include temperature as a variable in subsequent screens and optimization experiments.

Retain and observe plates until the drops are dried out. Crystal growth can occur within 15 minutes or one year.

Grid Screen MPD Formulation

Grid Screen MPD reagents are formulated using the highest purity chemicals, ultrapure water (18.2 Megohm-cm, 5 ppb TOC) and are sterile filtered using 0.22 micron filters into sterile containers (no preservatives added).

Grid Screen MPD reagents are stable at room temperature and are best if used within 12 months of receipt. To enhance reagent stability it is strongly recommended that Grid Screen MPD be stored at 4°C or -20°C. Avoid ultraviolet light to preserve reagent stability.

References and Readings

1. Crystallization of nucleic acids and proteins, Edited by A. Ducruix and R. Giege, The Practical Approach Series, Oxford Univ. Press, 1992.

2. Current approaches to macromolecular crystallization. McPherson, A. Eur. J. Biochem. 189, 1-23, 1990.

3. Protein and Nucleic Acid Crystallization. Methods, A Companion to Methods in Enzymology, Academic Press, Volume 1, Number 1, August 1990.

4. Advance in Protein Chemistry Volume 41. Pages 1-33 (Patricia C. Weber). Academic Press, 1991.

5. Current approaches to macromolecular crystallization., McPherson, A., Eur. J. Biochem. 189, 1-23, 1990.

6. Crystallization of Membrane Proteins. Edited by Hartmut Michel, 1990. CRC Press.

Technical Support

Inquiries regarding Grid Screen MPD reagent formulation, interpretation of screen results, optimization strategies and general inquiries regarding crystallization are welcome. Please e-mail, fax, or telephone your request to Hampton Research. Fax and e-mail Technical Support are available 24 hours a day. Telephone technical support is available 8:00 a.m. to 4:30 p.m. USA Pacific Standard Time.

> Hampton Research 34 Journey Aliso Viejo, CA 92656-3317 U.S.A. Tel: (949) 425-1321 • Fax: (949) 425-1611 Technical Support e-mail: tech@hrmail.com Website: www.hamptonresearch.com

© 1991-2018 Hampton Research Corp. all rights reserved Printed in the United States of America. This guide or parts thereof may not be reproduced in any form without the written permission of the publishers.



HR2-215 (pg 2)

								Tube	e (+/-)	-2-Methyl-2,4-pent	anediol T	ube	Buffer		
								#		[% v/v]		#			
				p	H										
								A1.	10		A	1. 0.1	M Citric acid	d pH 4.0	
		4	5	6	7	8	9	B1.	20		E	1. 0.1	M Citric acid	d pH 4.0	
	10	۸.1	10	12	Λ.Λ	٨٥	٨٩	C1.	40		C	1. 0.1	M Citric acid	d pH 4.0	
	10	AI	AZ	AJ	A4	AS	AO	D1.	65		E	1. 0.1	M Citric acid	d pH 4.0	
	00	D1	DO	Do	D/	DE	DC	Δ2	10		Δ	2 0 1	M Sodium a	cetate tribyd	rate nH 5.0
%	20	DI	DZ	БЗ	D4	DO	DO	R2	20		,	2 0 1	M Sodium a	cotato tribud	rate pH 5.0
<u> </u>	40	<u>C1</u>	02	C 2	C1	C 5	6	D2.	20			ו.0.1 ים חי	M Sodium a	cetate tribud	rate pH 5.0
MP I	40	U1	62	63	64	60	60	02.	40			$^{\prime 2.0.1}$	M Codium a	icelale trinyo	
	6E	D1	20	20	ПЛ	D5	DG	D2.	60		L	2. 0.1	M Sodium a	icelale trinyo	rate pH 5.0
	00	וט	DZ	03	04	05	00	A3.	10		A	3. 0.1	M MES mor	nohydrate pH	6.0
								B3.	20		E	3. 0.1	M MES mor	nohydrate pH	6.0
								C3.	40		C	3. 0.1	M MES mor	nohydrate pH	16.0
Tho	nЦ inc	licatod	00 00	oh Gric	I Saraa	n roog	ant ic	D3.	65		C	3. 0.1	M MES mor	nohydrate pH	l 6.0
i iie		licaleu	onea	ch Gho		n reay		A4.	10		A	4. 0.1	M HEPES p	H 7.0	
the <i>i</i>	ACTUA	NL рН	of the I	reagen	t at 25.	0°C. A	Шрн	B4.	20		E	4. 0.1	M HEPES n	H 7.0	
adju	stment	is have	e been	made	using I	Hydroc	nloric	C1	10		-	2/ 01	MHEPES		
acid	or Soc	dium h	ydroxid	e.				04. D4	65		с г	1. 0.1	MUEDES		
								D4.	10		L	- 0.1	MTRALLO	0	
								A5.	10		4	15. U.I	M INS PH 8.	0	
								В5.	20		E	5. 0.1	M Iris pH 8.	0	
								C5.	40		C	5. 0.1	M Tris pH 8.	.0	
								D5.	65		C	05. 0.1	M Tris pH 8.	.0	
								A6.	10		A	6. 0.1	M BICINE p	H 9.0	
								B6	20		F	6 0 1	M BICINE n	H 9 0	
								C6	40		Ċ	6 0 1	M BICINE p	HOO	
								D6	-0 65					U 0 0	
								D0.	60		L	0. 0.	IN DICINE P	9.0	
Ohamia		. .						tion Decoud							
Cnemic	ai An	aiysis	s and I	Recon	nmeno		otimiza	tion Reagent	S						
	7 10	00/ /./) O Ma		1	un a ali a l			NA 44						400.0
<u>HR2-62</u>	<u>27</u> - 10	0% (+/	-)-2-IVIE	etnyi-2,	4-penta	anedioi	(IMPD),	200 milliliters	Wr I	18.18 C ₆ H ₁₄ O	2 CAS NUI	nber [107-41-5]	EC NO 203-	-489-0
HB2-83	8 1 - 1 (M Cit	ric acid	l 100 r	nilliliter	s		Mr 19	2 13	$C_{e}H_{0}O_{7}$	CAS Nu	nber [77-92-91	EC No 201-	-069-1
<u></u>	<u>//</u> 1.0			, 1001		0			2.10	001007	0/10/110		// 02 0]	20110201	000 1
HR2-56	<u>69</u> - 1.0	M So	dium a	cetate	trihydra	ite, 100	millilite	rs M _r 136	6.08	CH ₃ COONa · 3I	H ₂ O CAS	Numl	per [6131-90-	-4] EC No	204-823-8
HP2-59	2 7 - 0 5		S mon	ohydra	to 100	millilite	ore	M. 21	3 25		240 04	Numh	or [1/522/-0	14-81 EC No	224-632-3
<u>11nz-30</u>	<u>m</u> - 0.0		.0 11011	lonyura	100		15		0.20	0611310040	120 040	Numb	61 [143224-3		224-032-3
HR2-58	35 - 1.0	MHE	PES, 1	00 mill	iliters			Mr 238	3.31	C8H18N2O4S	CAS Num	ber [7	365-45-9]	EC No 230-	907-9
										0 10 2 4					
HR2-58	<u>89</u> - 1.0	M Tris	s, 100 r	nilliliter	S			M _r 12 ⁻	1.14	C ₄ H ₁₁ NO ₃	CAS Num	ber [7	7-86-1]	EC No 201	-064-4
HP2-F0	10 - 1 0	MRIC		00 mill	ilitore			M. 160	3 17		CAS Num	har [1]	50-25-/1	EC No 205	-755-1
<u>nnz-30</u>	1 <u>3</u> - 1.0		JINE, I		111615			IVIT TOC	J. /	061131104		nei[]	0-20-4]	EC NU 205	-700-1

34 Journey Aliso Viejo, CA 92656-3317 U.S.A. Tel: (949) 425-1321 • Fax: (949) 425-1611 e-mail: tech@hrmail.com Website: www.hamptonresearch.com



Solutions for Crystal Growth

© 1991 - 2018 Hampton Research Corp. all rights reserved Printed in the United States of America. This guide or parts thereof may not be reproduced in any form without the written permission of the publishers.

Sample:	Sample Concentration:
Sample Buffer:	Date:
Reservoir Volume:	Temperature:
Drop Volume: Total μl Sample μl	Reservoirµl Additiveµl

_

	~	-	
1	Clear	Drop	

Microcrystals

2 Phase Separation 4 Birefringent Precipitate or

6 Needles (1D Growth) 3 Regular Granular Precipitate

7 Plates (2D Growth)

5 Posettes or Spherulites

8 Single Crystals (3D Growth < 0.2 mm)

9 Single Crystals (3D Growth > 0.2 mm)

Gı	id Screen [™] MPD - HR2-215 Scoring Sheet	Date:	Date:	Date:	Date:
A1.	0.1 M Citric acid pH 4.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol		1		
A2.	0.1 M Sodium acetate trihydrate pH 5.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol				
A3.	0.1 M MES monohydrate pH 6.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol	İ			
A4.	0.1 M HEPES pH 7.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol	1			
A5.	0.1 M Tris pH 8.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol				
A6.	0.1 M BICINE pH 9.0, 10% v/v (+/-)-2-Methyl-2,4-pentanediol				
B1.	0.1 M Citric acid pH 4.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
B2.	0.1 M Sodium acetate trihydrate pH 5.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
B3.	0.1 M MES monohydrate pH 6.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
B4.	0.1 M HEPES pH 7.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
B5.	0.1 M Tris pH 8.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
B6.	0.1 M BICINE pH 9.0, 20% v/v (+/-)-2-Methyl-2,4-pentanediol				
C1.	0.1 M Citric acid pH 4.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
C2.	0.1 M Sodium acetate trihydrate pH 5.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
C3.	0.1 M MES monohydrate pH 6.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
C4.	0.1 M HEPES pH 7.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
C5.	0.1 M Tris pH 8.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
C6.	0.1 M BICINE pH 9.0, 40% v/v (+/-)-2-Methyl-2,4-pentanediol				
D1.	0.1 M Citric acid pH 4.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				
D2.	0.1 M Sodium acetate trihydrate pH 5.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				
D3.	0.1 M MES monohydrate pH 6.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				
D4.	0.1 M HEPES pH 7.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				
D5.	0.1 M Tris pH 8.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				
D6.	0.1 M BICINE pH 9.0, 65% v/v (+/-)-2-Methyl-2,4-pentanediol				



34 Journey Aliso Viejo, CA 92656-3317 U.S.A. Tel: (949) 425-1321 • Fax: (949) 425-1611 e-mail: tech@hrmail.com Website: www.hamptonresearch.com

© 1991-2018 Hampton Research Corp. all rights reserved Printed in the United States of America. This guide or parts thereof may not be reproduced in any form without the written permission of the publishers.