

Data Sheet

Strep-Tactin[®] Sepharose[®]

Cat. No.: 2-1201-002, 2-1201-010, 2-1201-025,

2-1201-100, 2-1201-500

Lot No.: 1201-

IBA Headquarters IBA GmbH

Rudolf-Wissell-Str. 28 37079 Goettingen Germany Tel. +49 (0) 551-5 06 72-0 Fax +49 (0) 551-5 06 72-181

IBA US Distribution Center

1328 Ashby Road Olivette, MO 63132 USA Tel. 1-877-IBA-GmbH (1-877-422-4624) Fax 1-888-531-6813

E-mail: info@iba-lifesciences.com http://www.iba-lifesciences.com

Last date of revision March 2012 Version 1201-8

Description	Immobilized streptavidin variant called <i>Strep</i> -Tactin (5 mg/ml resin) which has been especially optimized for the purification of <i>Strep</i> -tag II fusion proteins*. Important note: To allow an efficient <i>Strep</i> -tag/ <i>Strep</i> -Tactin binding we strongly recommend using column purification instead of batch applications for proteins fused to <i>Strep</i> -tag II. It is crucial that protein binding takes place on the column. Even a pre-incubation of resin and lysate prior to filling the resin into a column will lead to decreased protein yields. Batch purification should be performed using <i>One-STrEP</i> -tag only. Further, prolonged batch incubations increase the risk of proteolytic degradation of the target protein including cleavage of the tag.
Support	Sepharose 4FF, 4% agarose
Form	50 % suspension in 100 mM Tris/HCl pH 8.0, 1 mM EDTA, 150 mM NaCl.
Biotin binding activity	> 300 nmol/ml resin
Stability	6 months after shipping
Storage	4 °C
Shipment	Room temperature

* Voss, S. & Skerra, A. (1997) Mutagenesis of a flexible loop in streptavidin leads to higher affinity for the *Strep*-tag II peptide and improved performance in recombinant protein purification. *Protein Eng.* 10, 975-982.

For research use only

Important licensing information

This product is based on Strep-Tactin technology covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at http://www.iba-lifesciences.com/patents.html or upon inquiry at info@iba-lifesciences.com or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

Trademark information