

Data Sheet

Strep-Tactin[®] Superflow[®] **high capacity**

Cat. No.: 2-1208-002, 2-1208-010, 2-1208-025,
2-1208-100, 2-1208-500

Lot No.: 1208-

Description	High capacity Strep-Tactin resin for the purification of Strep-tag II and One-STREP-tag fusion proteins. <i>Strep-Tactin</i> is a streptavidin variant with optimized binding properties for <i>Strep-tag</i> fusion proteins*. Important note: To allow an efficient <i>Strep-tag/Strep-Tactin</i> binding we strongly recommend using column purification instead of batch applications for proteins fused to <i>Strep-tag</i> II. It is crucial that protein binding takes place on the column. Even a pre-incubation of resin and lysate prior to filling the resin into a column will lead to decreased protein yields. Batch purification should be performed using <i>One-STREP-tag</i> only. Further, prolonged batch incubations increase the risk of proteolytic degradation of the target protein including cleavage of the tag.
Support	Superflow 6 (6% agarose, crosslinked)
Form	50 % suspension in 100 mM Tris/HCl pH 8.0, 1 mM EDTA, 150 mM NaCl
Biotin binding activity	> 900 nmol/ml resin
Elution	Buffer E (elution buffer): 100 mM Tris-Cl, 150 mM NaCl, 1 mM EDTA, 2.5 mM desthiobiotin, pH 8. It may be advantageous to use 5-10 mM desthiobiotin to get the target protein eluted at higher concentration.
Regeneration	HABA cannot be efficiently removed from Strep-Tactin Superflow High Capacity by using Buffer W. We recommend using Buffer W at pH 10.5 (or alternatively 100 mM Tris base) for efficient removal of HABA.
Stability	6 months after shipping
Storage	4 °C
Shipment	Room temperature

* Voss, S. & Skerra, A. (1997) Mutagenesis of a flexible loop in streptavidin leads to higher affinity for the *Strep-tag* II peptide and improved performance in recombinant protein purification. *Protein Eng.* 10, 975-982.

For research use only

Important licensing information

This product is based on Strep-Tactin technology covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at <http://www.iba-lifesciences.com/patents.html> or upon inquiry at info@iba-lifesciences.com or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

Trademark information

The owners of trademarks marked by "®" or "TM" are identified at <http://www.iba-lifesciences.com/patents.html>. Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.