



cellcrown







USER INSTRUCTIONS | MODE D'EMPLOI | GEBRAUCHSANWEISUNG | INSTRUCCIONES PARA EL USUARIO |  
ISTRUZIONI PER L'USO | BRUKSANVISNING |  
KÄYTTÖOHJE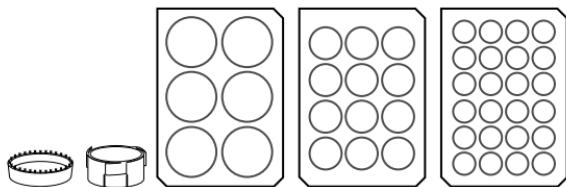

**A**

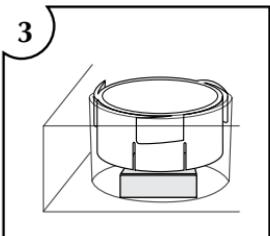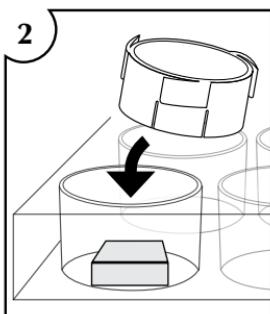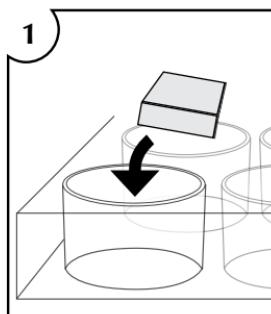**B**

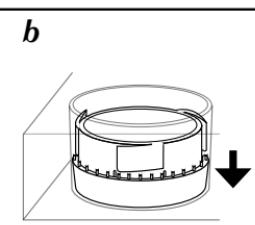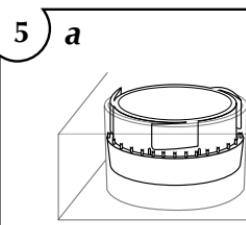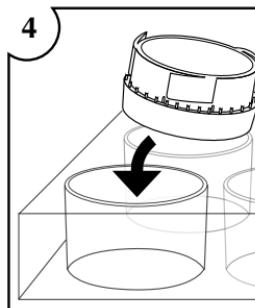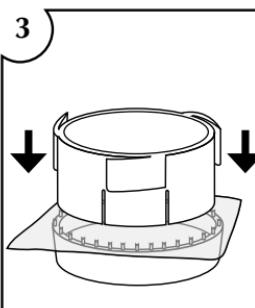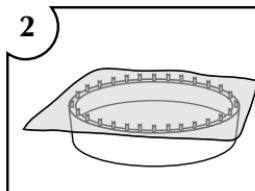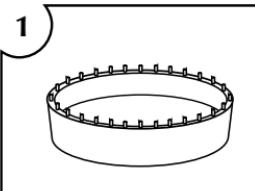EN	Sterilization temperature.....	121 °C / 250 °F
	Sterilization time.....	15 min
	Drying time.....	15 min

Table 1

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## CELLCROWN -INSERT

CellCrown is a disposable insert that immobilizes a sample in its place on the well plate. CellCrown makes it possible to combine various cell samples and materials in a cell culture environment, for example. The insert can be removed after culture.

CellCrown product selection includes sterile products ready for use and non-sterile products that will be sterilized by the customer. Inserts are manufactured for plates with 6, 12 and 24 wells (Table 2). CellCrown inserts fit to the most common standard well plates (Nunc, Corning). CellCrown inserts are made of polycarbonate (medical grade).

## USER INSTRUCTIONS

Figures A and B show ways of posing and using the CellCrown insert.

A

The CellCrown body can be used alone with rigid fibre, foil, block, textile, membrane and tissue samples, for example.

B

The combination of the CellCrown body and the ring can be used with flexible fibre, foil, textile, membrane and tissue samples. The height of the insert in the well is adjustable (Figure 5).

## STERILIZATION

CellCrown is disposable and must not be reutilized or re-sterilized. Scaffdex will not be responsible for any damage caused by any use contrary to instructions.

The CellCrown inserts that are sold non-sterile can be sterilized by gamma or beta radiation (with a maximum dose of 25 kGy), ethylene oxide gas, plasma or steam sterilization. The temperature and duration parameters tested and suitable for steam sterilization have been presented in table 1. Should the temperatures or durations indicated by the manufacturer be exceeded

for reasons of client's own sample, the insert might be damaged and become unusable.

Sterile CellCrown has been sterilized by gamma radiation (with a dose of 25 kGy).

A patent application for CellCrown has been registered.

Feedback and proposals are welcome! Our contact references are on the back page of the brochure.

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