



INTRODUCTION

Visikol® ENDO is a patent-pending chemical formula which quickly and easily renders animal tissues transparent. The Visikol® ENDO chemical formula is intended for use with large animal tissues and for gross visualization. Researchers have utilized Visikol® ENDO for many applications such as clearing whole mouse brains in conjunction with dye perfusion. Specific applications of Visikol® ENDO can be found on Visikol.com.

ADVANTAGES OF VISIKOL® ENDO

1. Rapid clearing time – 48 hours for whole mouse brain at 37°C.
2. Easy-to-use.
3. Non-destructive to underlying cellular structure of tissues.
4. No dehydration required.
5. Non-carcinogenic.

OPERATION

1. Visikol® ENDO is designed for use with whole mouse organs (e.g. mouse brain and kidney) or whole fetal animals (e.g. mouse, rat, rabbit).
2. A tissue can be fresh or fixed and is simply placed in Visikol® ENDO until clear.

Staining of tissues should take place prior to clearing and to accelerate the process tissues should be placed in a jar that is gently agitated and heated to 37°C.

Notice - Visikol® ENDO is not compatible with immunolabeling our fluorescent protein

STABILITY AND STORAGE

Visikol® ENDO can be stored at room temperature indefinitely, but should be kept out of direct lighting.

WARNINGS AND PRECAUTIONS

The MSDS for Visikol® ENDO can be found [here](#)

TECHNICAL ASSISTANCE

Email: order@visikol.com

MANUFACTURER

Visikol Inc.

www.visikol.com

120 Albany St, Tower II #850, New Brunswick, NJ 08901.

1-800-615-8474